

Cryostream Temperature Change Protocol

On the computer of the side you are working (left for the HF, left port and right computer for the HR, right port) open the program called ANAWIN (there is an icon on the desktop).

In the pull down menu called Channel Setup choose Channel Parameters. The Channel Setup screen will be displayed

Make sure the channel numbers selected is channel 1, if not use the arrow to select.

Highlight the number in the set point box and type in the value in centigrade you want the stream to be at and then hit enter.

The temperature should start to change as viewed in the process box below the set point.

You can also check the control box display below the detector.