


# Daily instrument cleaning guide

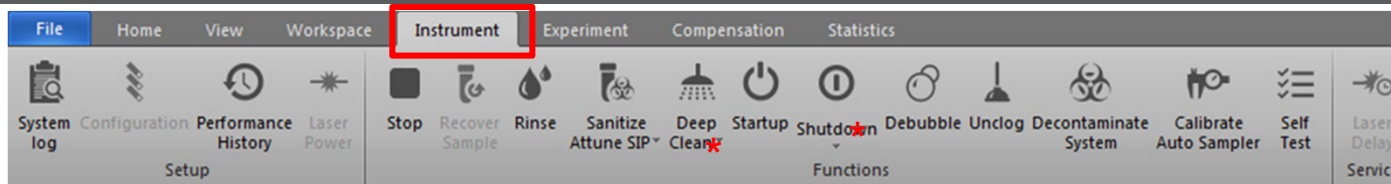

Between samples	<ul style="list-style-type: none"> <li>• <b>Rinse</b> – automatically initiated when SIP is lowered (for tubes), or set in <i>run protocol</i> for plates</li> <li>• <b>Sanitize SIP</b> between sticky samples or cell counts</li> </ul>
Between users / experiments	<ul style="list-style-type: none"> <li>• <b>Quick Deep Clean</b> - 30 minute cleaning routine (click on the arrow below the Deep Clean icon to select <u>Quick</u>)</li> <li>• <b>2x Sanitize SIP / Sanitize Autosampler SIP</b> (plate experiments) – <ul style="list-style-type: none"> <li>1<sup>st</sup> with 10% bleach per instructions</li> <li>2<sup>nd</sup> with MilliQ water.</li> </ul> </li> </ul>
End of day	<ul style="list-style-type: none"> <li>• <b>Thorough Shutdown</b> (click on the arrow below the Shutdown icon to select <i>Thorough</i>)</li> </ul>



*Note: Always wipe the outside of the SIP after doing a SIP Sanitize*